


Minifors 2


The Minifors 2 is a compact and easy-to-use bioreactor with a full range of application possibilities. It is a complete package that enables both, beginners and experienced users to easily perform applications. The Minifors 2 is available in two versions, one dedicated for microorganisms (M) and one dedicated for cell cultures (C).



Dimensions

Dimensions Overall Device

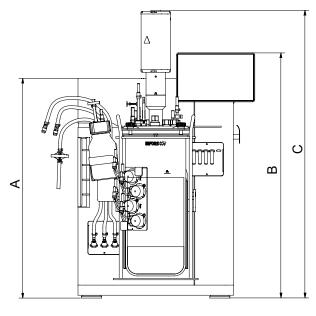


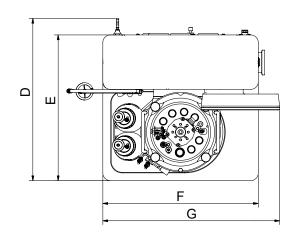



		Microorganisms (M)	Cell cultures (C)
А	Height basic unit without operating panel	565 mm	565 mm
В	Height basic unit incl. operating panel	631 mm	631 mm
С	Height incl. motor, vessels DN 145	740 mm	815 mm
	Height incl. motor, vessels DN 115	740 mm	815 mm
	Height incl. motor, vessels DN 90	770 mm	815 mm
D	Depth basic unit including hose nipples	415 mm	415 mm
Е	Depth basic unit without hose nipples	375 mm	375 mm
F	Width basic unit without operating panel	400 mm	400 mm
G	Width basic unit incl. operating panel	455 mm	455 mm



Dimensions Culture Vessel

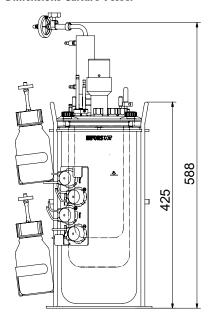

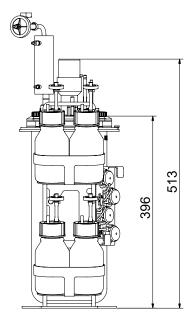

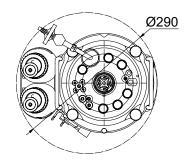

Weight

Basic unit	23.5 ± 0.5 kg

Culture Vessel

General	
Operating pressure in culture vessel	Pressureless
Form	Cylindrical with flat bottom
Material glass vessel	Borosilicate glass
Material top plate and built-in parts	AISI 316L, electropolished ¹⁾
Material O-rings (in contact with product)	EPDM

 $^{^{\}rm 1)}$ Exception: impellers in culture vessel 1.5 L / DN 90 for microorganisms are made of PEEK.

Vessel sizes			
Total volume [L]	1.5	3.0	6.0
Max. working volume [L]	1.0	2.0	4.0
Min. working volume version M [L]	0.3	0.6	1.1
Min. working volume version C [L]	0.3	0.7	1.5
Nominal diameter (DN, inner diameter vessel) [mm]	90	115	145
Height [mm]	235	295	370
Weight [kg] ¹⁾	6 ± 0.5	7 ± 0.5	9 ± 0.5

¹⁾ Equipped culture vessel, without medium, with vessel holder. The actual weight depends on design and allocation.

Ports in top plate		Quantity acc. to vessel DN		
Diameter	Thread	DN 90	DN 115	DN 145
7.5 mm	None	4	4	4
10 mm	None	4	4	4
12 mm	Pg13.5	4	6	7



Stirrer

General		М	С
Drive		Shaft with mechanical seal	
Direction of rotation of stirrer shaft		Counter-clockwise (top view)	
Bearing		Outside vesse	el, in drive hub
Motore type		DC, bri	ushless
Nominal power of	DN 90	102 W	74 W
motor	DN 115	260 W	74 W
	DN 145	260 W	74 W
Min. rotation speed		150 min ⁻¹	24 min ⁻¹
Max. rotation speed		1600 min ⁻¹	600 min ⁻¹
Accuracy	at ≤ 500 min ⁻¹	± 5 min ⁻¹	
measurement	at > 500 min ⁻¹	1 % setpoint	
Accuracy control		1% Full Scale	
Accuracy	at ≤ 300 min ⁻¹		± 2 min ⁻¹
measurement and control	at > 300 min ⁻¹		± 4 min ⁻¹

Impellers for microorganisms				
Туре		Rushton	impellers,	6 blades
Material impellers DN 145 and DN 115		AISI 316L, electropolished		
Material impellers DN 90		PEEK		
Quantity		2		
Dimensions impellers for microorganisms		DN 90	DN 115	DN 145
B		38 mm	46 mm	54 mm
	В	9 mm	11 mm	11 mm
ØA OA		11 mm	11 mm	11 mm

Impellers for cell cultures	
Туре	Pitched blade impeller with 3 blades, angled 45°
Material	AISI 316L, electropolished
Quantity	Standard: 1, option: 2
Flow direction blades	Standard: downwards, option: upwards

nm
nm
nm
n

Temperature Control System

Heating	Electrical, thermal block 630 W
Cooling	Tap water ¹⁾ via thermal block and adapter
Sensor	Type: Pt100 1/3 DIN-B
Measurement range	0 °C to +145 °C
Control range	Min. flow temperature + 5 °C to 60 °C
Accuracy measurement 2)	± 0.1 °C
Accuracy control 2)	± 0.2 °C

¹⁾ A circulating cooler can be used instead of tap water.

Gassing System

General specifications	М	С
Gas entry	Sparger	Sparger or head space (air and/or CO ₂ possible)
Specific gas flow rate 1)	20 L min ⁻¹	2000 mL min ⁻¹
Gas(ses)	Air, Air + O ₂ , Air + N ₂	Air, O ₂ , N ₂ , CO ₂
Flow control, mass flow controllers	2 pieces	5 pieces
Accuracy mass flow controllers	± 0.05 L min ⁻¹	± 4 mL min ⁻¹

¹⁾ Calculated for the max. working volume for all vessel sizes.

Control ranges of gas flow	М	С
Vessel 1.5 L TV / DN 90	0.05 to 2.0 L min ⁻¹	1.5 to 150 mL min ⁻¹
Vessel 3.0 LTV / DN 115	0.05 to 4.0 L min ⁻¹	3.0 to 300 mL min ⁻¹
Vessel 6.0 LTV / DN 145	0.05 to 8.0 L min ⁻¹	6.0 to 600 mL min ⁻¹

²⁾ +20 °C to +60 °C



pH Control

Control	
Control via cascade	Addition of acid and base via peristaltic pumps
	Version for cell cultures only: Addition of CO ₂ instead of acid possible
Control range	pH 2 to 12

Measurement system HAMILTON		
Sensor type Easyferm Plus ARC		
Measurement range pH 0 to 14		
Measurement system METTLER		
Sensor type	InPro3253i	
Measurement range	pH 0 to 12	

Pumps

General			
Туре		Peristaltic	
Quantity		4 pieces	
Control (operating modes)	Analogue	Continuous operation with variable speed	
	Digital	OFF/ON operation with fixed speed	

Hoses	Standard	Option 1	Option 2
Inside diameter	1.0 mm	0.5 mm	2.5 mm
Wall thickness	1.1 mm	1.15 mm	1.0 mm
Delivery rate 1)	3.5 ml min ⁻¹	1.1 ml min ⁻¹	16.1 ml min ⁻¹
Material	PharMed BPT		

¹⁾ Typical figure with water measured at max. rotation speed.

pO₂ Control

Control		
Control via cascade	Stirrer, gas flow, gas mixture (addition of O ₂ or N ₂)	
Control range	0 to 150 % O ₂ saturation	

Measurement system HAMILTON	
Sensor type	Visiferm DO ARC
Measurement range	0 % to 300 % air saturation

Measurement system METTLER	
Sensor type	InPro6860i, ISM
Measurement range	0 to 60 % O ₂ saturation

Operating Panel

НМІ	7" colour touch screen	
Operating system	Embedded Linux	
OPC server	OPC UA	

Antifoam Control

Sensor	Conductive with dosing needle
Control	Peristaltic pump <i>Antifoam</i>
Range	0 or 100 % (OFF or ON)

Turbidity Measurement (Optional)

Variant OPTEK		
Sensor type	ASD12-N	
Optical path lengths	OPL05 (higher cell densities) OPL10 (lower cell densities)	
Measurement range absorption	0 to 4 CU	

Variant aquila biolabs		
Sensor type	CGQ BioR	
Measurement modes	Green (521 nm) (low cell densities), Infrared (940 nm) (high cell densites)	
Measurement range	0 to 1000	



Exit Gas Analysis (Optional)

For aerobic bioprocesses	Measurement range	
	Vol. % O ₂	Vol. % CO ₂
Blue in One Ferm or Blue Vary	1.0 to 50	0 to 10
Blue in One Ferm or Blue Vary	1.0 to 50	0 to 25

For aerobic and anaerobic bioprocesses	Measurement range	
	Vol. % O ₂	Vol. % CO ₂
Blue in One Cell or Blue Vary	0 to 100	0 to 10
Blue in One Cell or Blue Vary	0 to 100	0 to 25

Interfaces

USB	USB 2.0 Type A Jack
Analog I/O	2x Input 4 to 20 mA 2x Output 4 to 20 mA
LAN	RJ45
Balance input 1)	9-pin D-SUB, RS232
Service interface	9-pin D-SUB, RS232

¹⁾ Supported balances (configuration by INFORS HT required): Kern DS 30K0.1, Kern FKB 6K0.02, Mettler MS32001L/01, Mettler MS6002TSDR/00

Redox Measurement (Optional)

Sensor type	Easyferm Plus ORP ARC
Measurement range	-1500 to +1500 mV

Various

IP rating	IP22
Sound pressure	< 70 dB (A)

Operating Conditions

Ambient temperature	5 °C to 40 °C
Ambient humidity	20 % to 90 %
Altitude operating location	max. 2000 m.a.s.l
Pollution degree as per EN 61010-1	2
Minimum distance	150 mm


Electrical Connection and Power Values

Mains voltage	120 / 230 VAC
Mains frequency	50 / 60 Hz
Max. power consumption 1)	~ 800 W
Max. current consumption	8 A
Fuse (two 5 x 20 mm fuses, time lag)	8 A

 $^{^{\}scriptsize 1)}$ During heating phase, vessel with max. 4 L working volume, at max. rotation speed.



Connections/Utilities

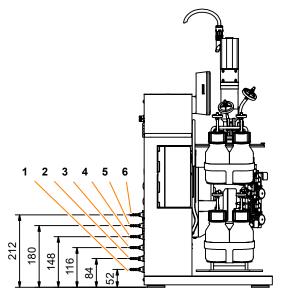

Pos.	Connection	Size	Pressure	Requirements
1	Water outlet	Hose nozzle 6 mm	No back pressure	
2	Water inlet	Hose nozzle 6 mm	2 ± 1 bar	■ Water hardness very soft or soft (CaCO ₃ concentration 0 to 1.5 mmol l ⁻¹) ■ Min. flow temperature: 10 °C
3	Air In	Hose nozzle 6 mm	2 ± 0.5 bar	■ Dry, clean and free of oil and dust
4	O ₂ /Gas2 In			Compressed air: Class 1,2,3,4 as per DIN ISO 8573-1
5	N ₂ In (C only)			
6	CO ₂ In (C only)			

eve®



eve® is a platform software for planning, execution and analysis of bioprocesses. eve® allows you to record bioprocess data and store it in a central database. The software offers workflows from simple bioprocesses to the planning and execution of complex strategies with various phases.

eve® makes it possible to generate and store bioprocess knowledge. Various libraries for storing information on organisms and culture media are available. Thanks to soft-sensors, additional knowledge can be generated.

In addition to INFORS HT products, biotech machines and analysis devices from third-part manufacturers can be connected. This makes it possible to holistically control, monitor and analyse bioprocesses using a single software.

eve* is installed on a centralised server. Access takes place via a browser, no client side installation is required. Bioprocess data is therefore available directly via the browser and independent of the operating system.

Various packages of the software are available. This makes it possible to adapt it to the individual needs and requirements of its users. eve® (in the premium version) is also suitable for working in a validated environment as per FDA CFR 21 Part 11.